

Effective Use of Various Aluminum Pans / Glass-Fiber Sheets

■ Aluminum pans

(1) AX-MX-30 Disposable Aluminum Pans



Ø 83 mm, 100 pcs/set

Suitable for samples that are sticky.

(2) AX-MX-31 Standard Aluminum Pans



Ø 85 mm, 100 pcs/set

Washable and reusable.

■ Glass-fiber sheets

(1) AX-MX-32-1 Glass-fiber Filtration Sheets



Ø 70 mm, 100 pcs/set

Suitable for liquid samples with high surface tension and high-sugar samples.

Can be used to cover dark samples so that they are not burnt.

* It is necessary to dry the sheet before use as it absorbs ambient moisture.

(2) AX-MX-32-2 Glass-fiber Sheets



Ø 78 mm, 100 pcs/set

Suitable for liquid samples, especially those that form a film/skin on the surface such as milk.

Can be used to cover dark samples so that they are not burnt.